


PAULA

Personal **AU**tomated **L**ab **A**ssistant

Release Notes 1.2.3

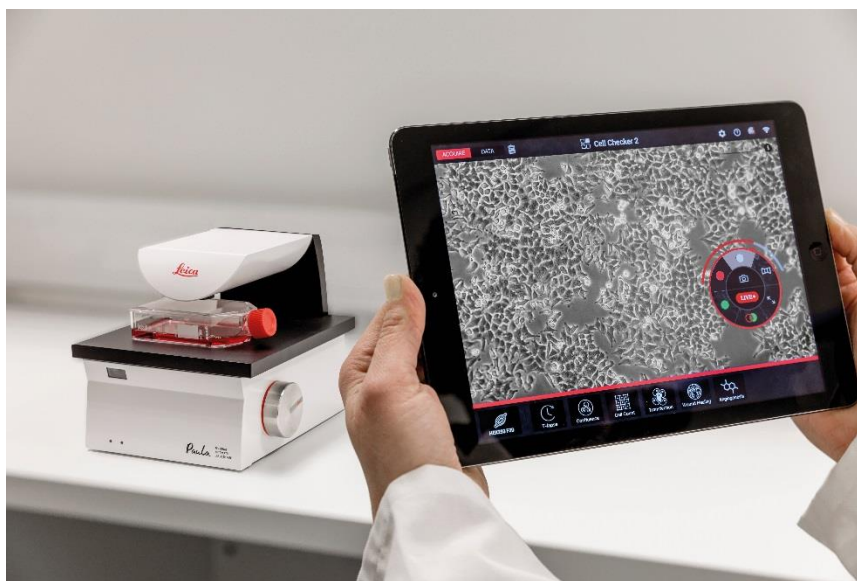

Release documentation for PAULA Patch 1.2.3

This document describes the 1.2.3 release of the PAULA imaging and analysis software for life science research. You should read this document before installing your copy of this software.

All reasonable steps have been taken to ensure that this publication is correct and complete, but should any user be in doubt about any detail, clarification may be sought from Leica Microsystems CMS GmbH, or their accredited representative. The information in this document is subject to change without notice and should not be construed, as a commitment by Leica Microsystems CMS GmbH.

Leica Microsystems CMS GmbH accepts no responsibility for any errors that may appear in this document.

Copyright © 2020 Leica Microsystems CMS GmbH

All rights reserved. The contents of this publication may not be reproduced in any form or communicated to a third party without prior written permission of Leica Microsystems CMS GmbH. Due to a policy of continuous development, we reserve the right to change specifications without notice. The term “Windows“, “Apple“, “iOS“, and “Android” may appear in the following text without further identification. The names of companies and products used herein may be trademarks of their respective owners.

Date: October, 2020

Content

1. Technical Requirements and released System Builds	4
1.1 Released System Builds.....	4
1.2 Minimum Requirements for Tablets and supported OS versions	4
2. Fixed Bugs/ Solved Issues	5
3. Recommendations	5
4. Open Issues/Restrictions.....	5

1. Technical Requirements and released System Builds

1.1 Released System Builds

Device	Released build
PAULA Service	1.2.3.26411
Android	1.2.2.26325
iOS	1.2.2.26318
UWP (Windows client)	1.2.2.26319
PAULA driver	1.2.3.14
PAULA firmware	V1.05
ELO Touch monitor driver	SW602623-EloMulti_Touch_6.9.17

Android and iOS can be updated via the stores if necessary. Windows PAULA client software (UWP) can be updated via PAULA webpage (Download tab).

1.2 Minimum Requirements for Tablets and supported OS versions

Device	Model	OS	Storage	RAM	Screen	Remarks
Android Tablet	Galaxy Tab S2	Support Android 8.0 and above	16 GB	3 GB	9.7-inch (1536 x 2048)	Minimum Android Version 8.0 required for downloading from Play Store and Baidu .
	Galaxy Tab A (2016) with S Pen					
	Galaxy Tab S3					
	Galaxy Tab S4					
iPAD	iPad Air 2	IOS Version 12.1.4 and above	32 GB	2 GB	9.7-inch (1536 x 2048)	
	iPad 6 th Gen					
	iPad Pro					
Window Device	Microsoft Surface	Windows 10 Pro Version 1803 (OS Build 17134.407)		8 GB	10.6-inch (1366 x 768)	Processor: Intel Core i7-6650u CPU@ 2.20GHz 64 bit OS x64-based processor, Window Defender Firewall - Enable

2. Fixed Bugs/ Solved Issues

- Update of license components based on WIBU Systems to a new version

3. Recommendations

- For an optimal performance we strongly recommend to install the released builds and to update your client (tablet or windows client of the NUC) if necessary with the recommended versions.
- We recommend to insert a valid e-mail address so you can retrieve a new password when your current password is forgotten or you typed 5 times a wrong password.

4. Open Issues/ Restrictions

- “Remote Care” cannot inform users proactively about issues
- To view TIFF formatted images you may need to install a 3rd party app on your tablet